

Cutting bands directly out of UV excited gels

We do not allow this option due to a number of issues:

1. Health and Safety concerns
2. Damage to gel tray/box




HOWEVER we have a far better approach to negate these two issues...

Amersham Imager Solution

1. Image gel
2. Print out 1:1 image using software
3. Lay gel on top of image (either print to cellulose acetate or place glass/plastic support between gel and paper)
4. Cut out bands
5. Re-image gel to check bands excised correctly

Now can avoid “friendly lab. messages”

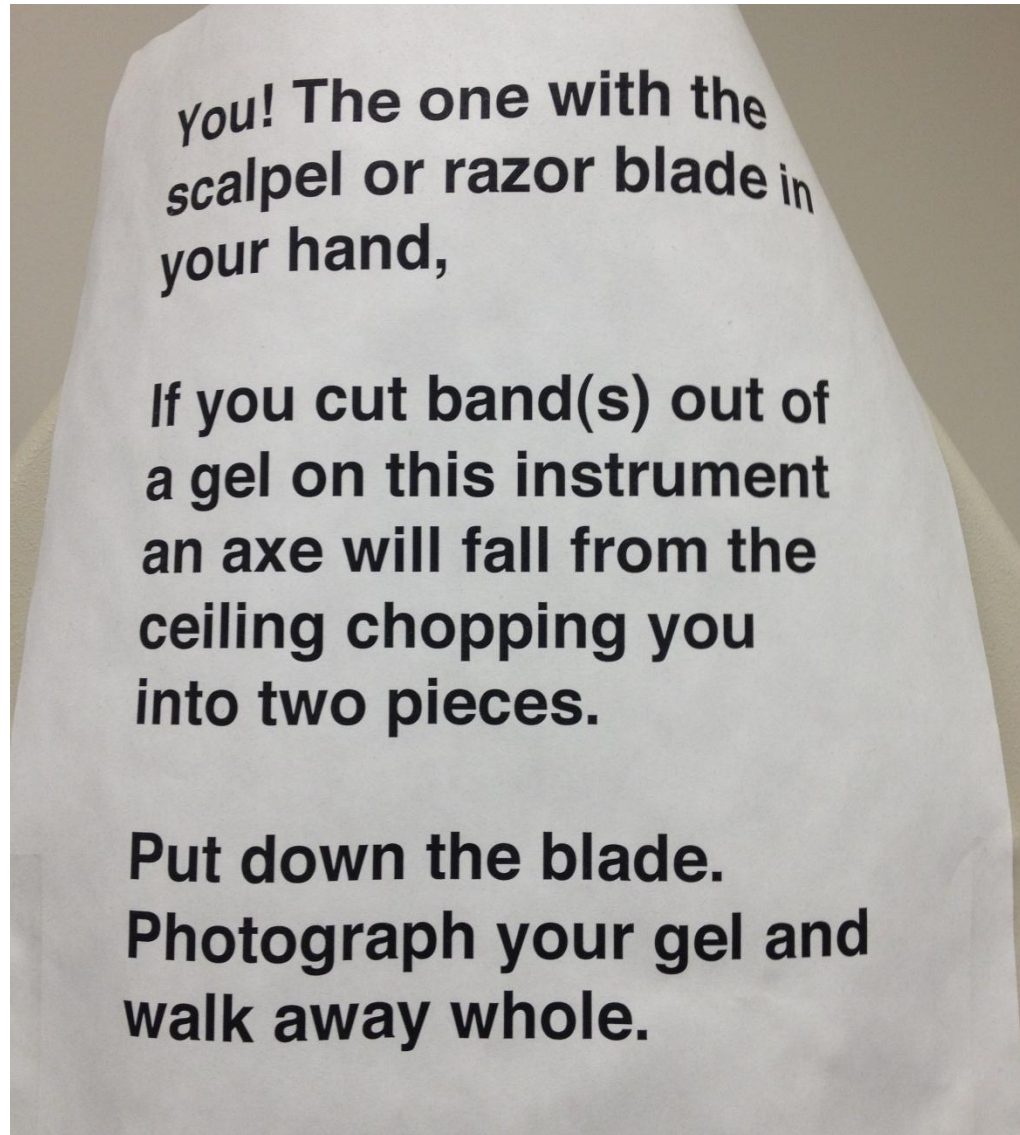



imagination at work